

Grounding of the Axioskop FS

Installation Instructions

Attach grounding connector (11) to the microscope:

Use Philips screw of the lower cover plate on the back of the microscope. The fixation point (2) should preferably be chosen on the bottom right.

Bonding of the lower part of the stand with the upper part:

Attach short cable (9) to the cover plate at (1) with Philips screw (10). Insert sprocket washer (8).

Bonding of the condenser carrier with the stand

Attach connecting piece (5) with washer (6) for the screwhead and sprocket washer (4) for the condenser carrier to the rear drilled hole (13) of the condenser carrier from below. Slide plug-on sleeve of the medium-long cable on the connecting piece.

If the microscope is equipped with a polarizer carrier, the stop bolt can be used for attachment of the connecting piece (5) instead of the screw (6). Washer (4) must be inserted in any case.

Grounding to the microscope:

Attach long cable and cable terminal (3) to an external grounding point with a screw.

Note:

Securely tighten all screws to make the microscope potential-free.

